



Smart, safe and clean chemistry

Gas Module

User's Manual

Version: 2.00

Date: June, 2013

1. Exchange the mixer

The mixer can be replaced by removing the top plate (see Figure 1.1). Next remove the mixer by disconnecting the mixer as depicted in Figure 1.2.



Figure 1.1: Disconnection of the top plate.

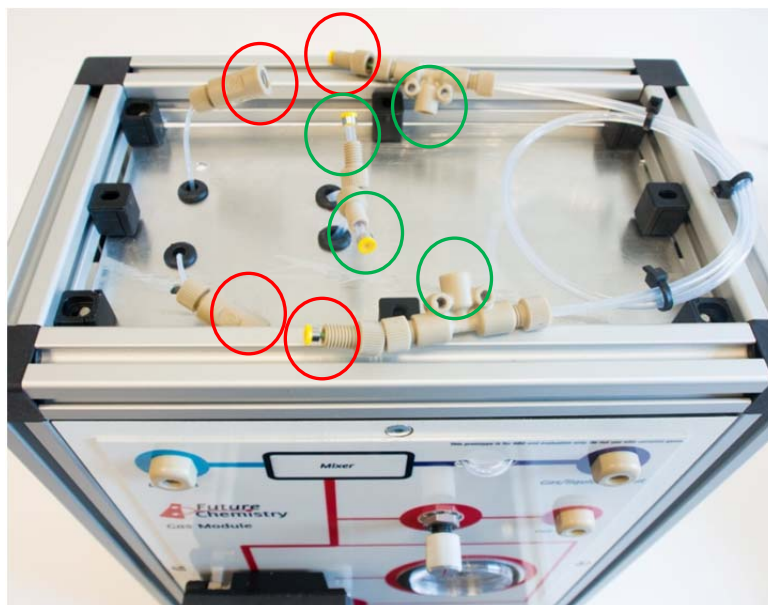


Figure 1.2: Disconnection of the mixer.

- WARNING -



CAUTION: Disconnect the mixer **ONLY** at the positions as depicted in Figure 1.2



FutureChemistry Holding BV
Toernooiveld 100
6525 EC Nijmegen
The Netherlands
Tel +31 24 711 4029
Fax +31 24 89 1446
E-mail: support@futurechemistry.com

© COPYRIGHT BY FUTURECHEMSITRY HOLDING BV